VFC2000-MT

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Device Maintenance

Battery Replacement

Materials: U9VL-J Battery or any 9 V Battery (lithium recommended)

1. On the bottom of the data logger, open the battery compartment by pulling in on the cover tab.
2. Remove the battery by pulling it from the compartment.
3. Install the new battery, taking note of the polarity.
4. Push the cover closed until it clicks.

Recalibration

Recalibration is recommended annually or bi-annually for any data logger; a reminder is automatically displayed in the software when the device is due. To send devices back for calibration, visit **[madgetech.com](https://www.madgetech.com)**.

Product Support & Troubleshooting:

- Refer to the Troubleshooting section of this document.
- Visit our Knowledge Base online at **[madgetech.com/resources](https://www.madgetech.com/resources)**.

EO: sales@eofirm.com | www.eofirm.com | 0755-84200058

MadgeTech 4 Software Support:

- Refer to the built-in help section of the MadgeTech 4 Software.
- Download the MadgeTech 4 Software Manual at **[madgetech.com](https://www.madgetech.com)**.

EO: sales@eofirm.com | www.eofirm.com | 0755-84200058

PRODUCT USER GUIDE

Specifications

EO: sales@eofirm.com | www.eofirm.com | 0755-84200058

TEMPERATURE	
Temperature Range	-20 °C to +60 °C (-4 °F to +140 °F)
Resolution	0.01 °C (0.018 °F)
Calibrated Accuracy	±0.50 °C/± 0.18 °F (0 °C to +55 °C/32 °F to 131 °F)
Response Time	10 minutes free air

REMOTE CHANNEL	
Thermocouple Connection	Female subminiature (SMP) (MP model) Pluggable screw terminal (TB model)
Cold Junction Compensation	Automatic, based on internal channel
Max. Thermocouple Resistance	100 Ω
Thermocouple K	Included Probe Range: -100 °C to +80 °C (-148 °F to +176 °F) Glycol Bottle Range: -50 °C to +80 °C (-58 °F to +176 °F) Resolution: 0.1 °C Accuracy: ±0.5 °C
Response Time	τ = 2 minutes to 63% of change

GENERAL	
Reading Rate	1 reading every second up to 1 reading every 24 hours
Memory	16,128 readings
LED Functionality	3 status LEDs
Wrap Around	Yes
Start Modes	Immediate and delay start
Calibration	Digital calibration through software

Calibration Date	Automatically recorded within device
Battery Type	9 V lithium battery included; user replaceable with any 9 V battery (lithium recommended)
Battery Life	3 years typical at 1 minute reading rate
Data Format	For Display: °C or °F For Software: Date and time stamped °C, K, °F or °R
Time Accuracy	± 1 minute/month
Computer Interface	USB to mini USB, 250,000 baud for standalone operation
Operating System Compatibility	Windows XP SP3 or later
Software Compatibility	Standard Software version 4.2.21.0 or later
Operating Environment	-20 °C to +60 °C (-4 °F to +140 °F), 0 %RH to 95 %RH non-condensing
Dimensions	3.0 in x 3.5 in x 0.95 in (76.2 mm x 88.9 mm x 24.1 mm) Data logger only
Glycol Bottle	30 mL
Probe Length	72 in
Material	ABS Plastic
Weight	4.5 oz (129 g)
Approvals	CE
Alarm	User configurable high and low audible and on-screen alarms. Alarm Delay: A cumulative alarm delay may be set in which the device will activate the alarm (via LED) only when the device has recorded a user specified time duration of data.
Audible Alarm Functionality	1 Beep per second for reading alarm above/below threshold

BATTERY WARNING: BATTERY MAY LEAK, FLAME OR EXPLODE IF DISASSEMBLED, SHORTED, CHARGED, CONNECTED TOGETHER, MIXED WITH USED OR OTHER BATTERIES, EXPOSED TO FIRE OR HIGH TEMPERATURE. DISCARD USED BATTERY PROMPTLY. KEEP OUT OF REACH OF CHILDREN.

Ordering Information

VFC2000-MT	PN 902311-00	VFC Temperature data logger with thermocouple probe and USB to mini USB cable
VFC2000-MT-GB	PN 902238-00	VFC Temperature data logger with thermocouple probe, glycol bottle and USB to mini USB cable
Power Adapter	PN 901839-00	Replacement USB universal power adapter
U9VL-J	PN 901804-00	Replacement battery for VFC2000-MT

■Tel:+86 755-8420 0058 ■Fax:+86 755-2822 5583 ■E-mail:sales@eofirm.com ■http://www.eofirm.com